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apotec[®] cool connect pharmaceutical refrigerator
in acc. with DIN 58345

USER MANUAL

apotec[®]

The professional solution for your lab.

COPYRIGHT

The copyright for this operating manual is owned by WEPA Apothekenbedarf GmbH & Co KG. This manual must not be reproduced in any format without the prior written permission of WEPA Apothekenbedarf GmbH & Co KG. WEPA Apothekenbedarf GmbH & Co KG shall not be liable for use of the information contained in these instructions. Errors excepted.

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TABLE OF CONTENTS

Foreword	4
Technical data	5-6
Safety information	7-12
Explanation of the terminology, operating states and alarm notifications	13
Possible error notifications	14
Warranty and liability	15
Notes	16
Warranty card	17

FOREWORD

The apotec® cool connect pharmaceutical refrigerator stores medicinal products requiring refrigeration and fulfils all requirements in accordance with DIN 58345, such as:

- Integrated temperature recording
- Safety thermostat
- Visual and acoustic alarm notification
- Battery for alarm in the case of power outage (12 h battery)

The device fulfils DIN 58345 only with the factory temperature setting of 5 °C. If a temperature adjustment occurs (can only be performed by the service technician) conformity for the purpose of DIN 58345 is no longer fulfilled.

At the same time, the apotec® cool connect pharmaceutical refrigerator in accordance with DIN 58345 from WEPA Apothekenbedarf offers you further product advantages, such as:

- Integrated data recording and overview on the spacious 7" display
- A potential-free relay contact for alarm relaying
- Convection cooling (see page 9)
- Automatic defrosting and evaporation of defrost water (see page 12)
- Simple menu navigation
- Alarm notifications possible by SMS or email

TECHNICAL DATA (Equipped with plastic drawers)

Cooling content:	236 l
Order no.:	019891 (door hinge, right), 019890 (door hinge, left)
Mains connection:	230 V
Size:	600 x 1365 x 610 mm (BxHxD)
Weight	65 kg
Drawers:	5 plastic drawers, max. load 15 kg per drawer (4 drawers with 2 lengthwise partitions and 6 crosswise partitions, 1 drawer with 1 lengthwise partition and 4 crosswise partitions)
Temperature display range:	-2 °C to +25 °C, maximum and minimum temperature display
Power consumption:	0.8 kWh / 24 h

Cooling content:	335 L
Order no.:	019894 (door hinge, right), 019893 (door hinge, left)
Mains connection:	230 V
Size:	600 x 1775 x 610 mm (BxHxD)
Weight	83 kg
Drawers:	9 plastic drawers, max. load 15 kg per drawer (8 drawers with 2 lengthwise partitions and 6 crosswise partitions, 1 drawer with 1 lengthwise partitions and 4 crosswise partitions)
Temperature display range:	-2 °C to +25 °C, maximum and minimum temperature display
Power consumption:	0.9 kWh / 24 h

Cooling content:	335 l – with glass door
Order no.:	019897 (door hinge, right), 019896 (door hinge, left)
Mains connection:	230 V
Size:	600 x 1775 x 610 mm (BxHxD)
Weight	90 kg
Drawers:	9 plastic drawers, max. load 15 kg per drawer (8 drawers with 2 lengthwise partitions and 6 crosswise partitions, 1 drawer with 1 lengthwise partition and 4 crosswise partitions)
Temperature display range:	-2 °C to +25 °C, maximum and minimum temperature display
Power consumption:	1,26 kWh / 24 h

Suppliable accessories up to 335 l: Order no.: 019867 crosswise partition 136 mm (VE 10 St.)
 Order no.: 019868 crosswise partition 206 mm (VE 10 pcs.)
 Order no.: 019869 card pockets (VE 10 St.)

TECHNICAL DATA (Equipped with AluCool drawers)

Cooling content: 236 l
Order no.: 019922 (door hinge, right), 019923 (door hinge, left)
Mains connection: 230 V
Size: 600 x 1365 x 610 mm (BxHxD)
Weight 65 kg
Drawers: 5 AluCool drawers, max. load 15 kg per drawer (all drawers with 2 lengthwise partitions and 6 crosswise partitions)
Temperature display range: -2 °C to +25 °C, maximum and minimum temperature display
Power consumption: 0.8 kWh / 24 h

Cooling content: 335 L
Order no.: 019920 (door hinge, right), 019921 (door hinge, left)
Mains connection: 230 V
Size: 600 x 1775 x 610 mm (BxHxD)
Weight 83 kg
Drawers: 9 AluCool drawers, max. load 15 kg per drawer (all drawers with 2 lengthwise partitions and 6 crosswise partitions)
Temperature display range: -2 °C to +25 °C, maximum and minimum temperature display
Power consumption: 0.9 kWh / 24 h

Cooling content: 335 l – with glass door
Order no.: 019924 (door hinge, right), 019925 (door hinge, left)
Mains connection: 230 V
Size: 600 x 1775 x 610 mm (BxHxD)
Weight 90 kg
Drawers: 9 AluCool drawers, max. load 15 kg per drawer (all drawers with 2 lengthwise partitions and 6 crosswise partitions)
Temperature display range: -2 °C to +25 °C, maximum and minimum temperature display
Power consumption: 1,26 kWh / 24 h

Suppliable accessories up to 335 l: Order no.: 600052 lengthwise partitions
Order no.: 600055 crosswise partitions
Order no.: 019869 card pockets (VE 10 St.)
Order no.: 600058 drawer for large packs
Order no.: 600059 lockable drawer

Cooling content: 572 l
Order no.: 019798 (door hinge, right), 019799 (door hinge, left)
Mains connection: 230 V
Size: 750 x 1775 x 730 mm (BxHxD)
Weight 92 kg
Drawers: 9 AluCool drawers, max. load 15 kg per drawer (all drawers with 3 lengthwise partitions and 8 crosswise partitions)
Temperature display range: -2 °C to +25 °C, maximum and minimum temperature display
Power consumption: 0,983 kWh / 24 h

Suppliable accessories 572 l: Order no.: 600053 lengthwise partitions
Order no.: 600055 crosswise partitions
Order no.: 019869 card pockets (VE 10 St.)
Order no.: 600056 drawer for large packs
Order no.: 600057 lockable drawer

SAFETY INFORMATION

Before using the refrigerator, read through these simple instructions. Non-compliance might have dangerous consequences or contravene regulations. The manufacturer shall not be liable for any damage arising from improper use of the equipment.

In general:

Check to see there is no damage to the refrigerator on unpacking it. Only operate the refrigerator as specified in the manual to ensure failsafe operation.

Disposal information:

The packaging is made of recyclable materials.

- Corrugated cardboard / cardboard
- Moulded parts made of foamed polystyrene
- Polyethylene film
- Polypropylene straps
- **Do not allow children to play with packaging material – films pose a risk of suffocation!**
- Take the packaging to an official collection point.

Disposal

The facility is made predominantly of steel (also of aluminium to a certain extent) (except for the electrical equipment) and must be disposed of according to the applicable local environmental regulations. Dispose of according to constitution, existing regulations and heeding current conditions, e.g. as:

- electrical waste (printed circuit boards);
- batteries, fluorescent lamps/energy-saving bulbs, hazardous substances;
- plastics (housing);
- sheet metal, steel, copper, aluminium (separated by metal type).

Before disposing of the device, decontaminate all parts that have been in contact with process substances. Remove any hazardous materials from the the apotec® pharmaceutical refrigerator.

Oils, solutions and detergents and contaminated cleaning tools (brushes, cloths, etc.) must be disposed of according to local conditions according to applicable waste codes and heeding the information found in the manufacturer's data sheets.

Battery regulation:

Do not dispose of the batteries in normal household waste. You are legally obliged to ensure that batteries are processed by a specialist disposal company. For this purpose, you can take your old batteries to the public collection points in your district free of charge or anywhere where batteries are sold. Alternatively, you can return the batteries to us at the following address

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We will then dispose of them properly and free of charge for you. Batteries containing toxic substances are marked with a symbol, consisting of a crossed out dustbin and the chemical symbol (Cd, Hg or Pb) of the heavy metal crucial for the toxic classification. The crossed out dustbin symbol indicates batteries must not be discarded in household waste. The chemical symbols indicate the following: Pb: battery contains lead; Cd: battery contains cadmium; Hg: battery contains mercury.

Place of installation / Power connection:

- Do not set up appliance in direct sunlight, near to heating and similar
- The floor must be level and horizontal at the location.
Correct uneven sections using adjustable feet.
- Ensure there is adequate ventilation and air circulation
- Do not use an extension lead to connect the refrigerator to other appliances – risk of over-heating!
- The place of installation must have a temperature range from 10 °C to 35 °C. The relative humidity may not exceed 75%.

Power connection:

Type of current (AC) and voltage at the point of installation must be the same as the particulars on the type plate. The type plate is inside the refrigerator on the left-hand side. The socket must be fused at 10 amperes or higher, be situated away from the rear side of the appliance and be easily accessible. The appliance must only be connected via a properly earthed earthing contact socket.

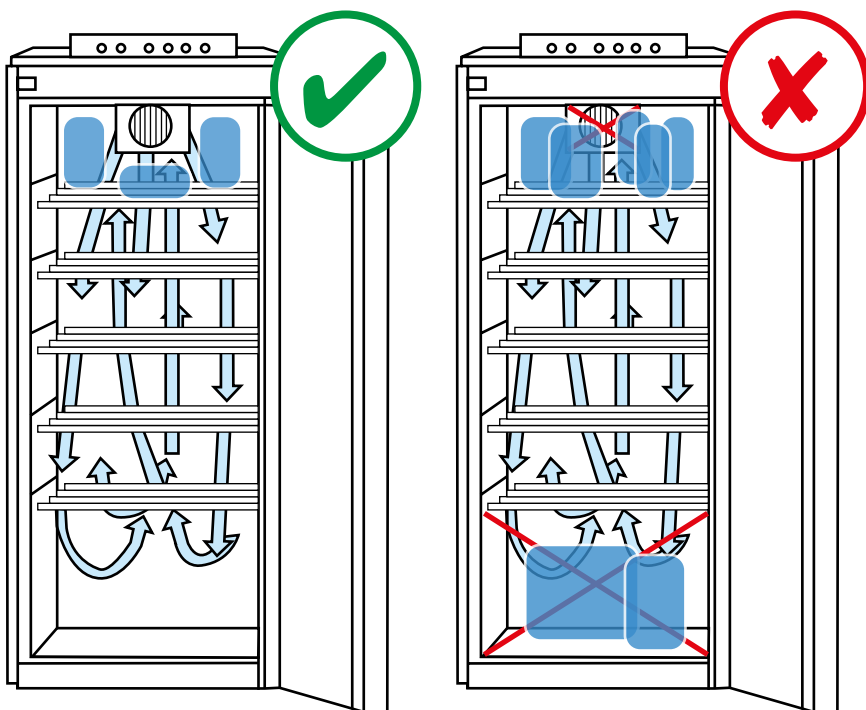
Recirculated air cooling:

Recirculated air cooling involves the swirling of cooled air with a fan. This process enables the cold air to be rapidly and evenly spread inside. Temperature stratification practically does not occur. The temperature is roughly equal throughout the interior.

Safety and warning information:

- To avoid any injuries or material damage, two people should be entrusted with unpacking and setting up the refrigerator.
- If any damage is noticed on the refrigerator, contact the supplier immediately – before connecting the appliance.
- To ensure failsafe operation, only install and connect the refrigerator in accordance with the instructions in the manual.
- If a fault occurs, disconnect the refrigerator from the mains, pull out the mains plug or trip/remove the fuse. **To disconnect the refrigerator from the mains, pull the mains plug itself and not the connecting cable.**
- **Interventions and repairs** to the refrigerator must be carried out by the competent Customer Services department. Failure to do so creates considerable risks for the user. The same applies to the replacement of the power connection line.
- Do not work with naked flames or ignition sources **inside the appliance**.
- Ensure that the refrigeration circuit is not damaged when transporting or cleaning the refrigerator. Keep sources of ignition away from any damaged points and ventilate the area well.
- Do not use **base, drawers, doors etc.** for stepping on or for propping up.
- This refrigerator is not intended for use by people (including children) with physical, sensory or mental impairments or for those who are inexperienced in its use or who lack the necessary knowledge – unless they have been instructed in using the refrigerator by someone who is responsible for their safety or who supervises them in the initial stages. Children should be supervised at all times to ensure they do not play with the refrigerator.
- Avoid constant skin contact with cold surfaces or cooled/frozen products. This can otherwise lead to pain, numbness and frost-bite. Provide for protection if work involves lengthy periods of skin contact e.g. by wearing gloves.

- Do not store in the refrigerator any explosive materials or spray cans with combustible propellants such as propane, butane and pentane. Any escaping gases could be ignited by electrical components. These kinds of spray cans can be recognized as such by the printed details on content or a flame symbol
- Do not use any electrical equipment inside the refrigerator.
- Keep keys of lockable equipment well out of the reach of children.
- Only load the refrigerator with 2 - 8°C refrigerated products (intrinsic temperature). Proper cooling cannot be ensured if this is disregarded.
- Do not put any foodstuffs/drinks into the refrigerator!
- Visually check information labels, type plate and magnetic contact at least twice a year as to their legibility/intactness. If any damage is detected, please immediately contact the Customer Services department.
- Do not keep any „frozen products“ (cold packs/pads, ice etc.) in the refrigerator. Proper cooling cannot be ensured if this is disregarded!
- To ensure a proper data record, date and time must always be up to date. Please check this regularly.

The lower/upper section of the refrigerator (see drawing) must not be blocked.
The recirculation function will be disturbed and proper ventilation cannot be guaranteed!

- The mains plug must always be freely accessible!
- The device is not suitable for installation, e.g. in a cabinet.
- The electronic unit with display must be easily accessible, to be able to operate the rear panel if necessary (connections, antenna, switches).

WARNING:

Risk of injury due to electric shock! The device may only be opened by the customer service or specially trained personnel.

- According to the EN 378 standard, the room where your device is installed must have a volume of 1 m³ per 8 g of refrigerant R 600a, so that no ignitable gas-air mixture can form in the installation room in the event of a leak in the refrigerant circuit. The refrigerant quantity is indicated on the rating plate inside the device.

CAUTION:

The device must be aligned horizontally and vertically. If the device is tilted, the body can become deformed and the door will not close properly.

- Level out uneven floors with the adjustable feet.
- Do not connect the device via extension cables or distribution sockets.

Fault:

You can yourself correct the following faults by checking on the possible causes:

- **If the appliance is not working:** Please check that
 - the appliance is switched on
 - the mains plug is plugged into the socket correctly,
 - the fuse on the plug is working.
- **If noises are too loud:** Please check that
 - the appliance is standing firmly on the floor,
 - the operating refrigerating unit is the reason for vibrations in furniture or objects next to it.Note that flow noises in the refrigeration circuit cannot be prevented.

If none of the aforementioned causes are present and you are unable to rectify the defect yourself, please contact the Customer Services department.

Defrosting:

The cooling section defrosts automatically.

If the rear wall still freezes over, the appliance must be defrosted manually.

- To defrost, switch off the refrigerator. Disconnect the mains plug
- Remove the cooling products from the refrigerator and store in a cool place.
- Leave the appliance door open when defrosting. Mop up any condensed water with a cloth and clean the refrigerator.

When defrosting, do not use any mechanical implements or other artificial aids save for those recommended by the manufacturer.

Cleaning:

- Switch the refrigerator off before cleaning (pull out the mains plug).
- Clean the interior, fitments (drawers) and external panels with tepid water and a little rinsing agent. Do not use any cleaning or chemical solvents containing sand or acid.
- Do not work with steam cleaning devices. Risk of damage/injury.
- Make sure that no cleaning water gets into the electrical parts or into the ventilation grate.
- The chiller with the heat exchanger (metal grate at the back of the refrigerator) should be cleaned or dusted once a year.
- Clean the drain in the refrigerator with something thin (Q-tips and the like).
- The operating display may only be cleaned with a soft, dry microfibre cloth. Under no circumstances use cleaning agents, alcohol, etc. Non-compliance can lead to irreparable damage to the display!

Battery:

- All rechargeable batteries wear out over time and have a limited life. The wear depends on many factors such as charging cycles, temperature influence, use, etc. We therefore recommend that the batteries be replaced by an authorised specialist company after 36 months at the latest.

Screensaver:

- Your device has a screensaver that starts after 2 minutes of inactive operation. As long as the screensaver is active, all functions are o.k. and the temperature is within the setpoint range (see also green LED). As soon as an error occurs, the screensaver automatically goes off. Otherwise, the screensaver can of course also be deactivated by touching the display.

EXPLANATION OF THE TERMINOLOGY, OPERATING STATES AND ALARM NOTIFICATIONS

Internal temperature:

Currently measured air temperature in the refrigerator.

Reference temperature:

„Simulates“ the temperature of the refrigerated products. This was recorded in accordance with DIN 58345 in a reference body facility. This temperature is also crucial for the alarm.

Target temperature:

The target temperature is preset in accordance with DIN 58345 at 5 °C. A deviation of ± 3 °C is permitted. The deviation refers to the reference temperature. If the temperature range is exceeded or fallen short of, the refrigerator emits a visual and acoustic alarm.

Note:

You will notice a short-term temperature increase every day which is caused by the automatic defrost function. However, the 8 °C reference temperature is not normally exceeded.

LEDs:

In addition to the indicated errors or warnings, the LED light shows an overview of the operating status:

Green:

Proper refrigerator operation

Yellow:

Appliance had an error but is in an operational state again (see error description in display). The error must be confirmed by the "Alarm Quitt" button. Please check the temperature profile /error report.

Red:

Appliance is not in an operational state and has a defect. See error notification in display. Steps must be taken without delay.

Start-up:

The refrigerator is in the non-operating mode at the time of delivery: the display is non-active. Inserting the mains plug deactivates the non-operating mode. The appliance incorporates a battery which has been charged for 36 hours on first-time operation.

Menu:

You can make the following settings in the menu.

- Personalisation
- Language selection
- Warning notifications
- Time setting
- Data export
- Temperature display
- Set up W-LAN connection
- Alarm volume
- Software
- Service (reserved for the Customer Services department)

POSSIBLE ERROR NOTIFICATIONS

Power failure:

If the appliance is disconnected from the mains cable or is no longer receiving power due to another cause (e.g. power outage) the **“power failure” error notification appears in the display and the background lighting is dimmed.** (Note: The display remains in this state for min. 12 hours or until the power supply is resumed). The green light, which indicates proper operation of the appliance, is extinguished. In the interim, the red LED starts to light up and the installed buzzer commences a signal. In this case, the buzzer can be switched off with the **“Alarm off”** button. When the power supply or mains voltage is switched on again, the error notification disappears after acknowledgement and normal operation is resumed. The device is in emergency mode and is switched back to normal mode by acknowledgement.

Caution: Only possible when the mains voltage is present again.

Exceeding and falling short of the reference temperature:

If the factory set 5 °C is exceeded or fallen short of by 3 °C, the red LED starts to light up and the buzzer sounds. The buzzer can be switched off with the **„Alarm off”** button. The **“reference temperature”** notification appears on the display. The notification and the red light remain active until the refrigerator re-enters the target range. The device is in emergency mode and is switched back to normal mode by acknowledgement.

Caution: Only possible when the device is cooling in the setpoint range again.

Door open:

If the door is open for longer than 120 seconds, an acoustic alarm sounds and an alarm notification also appears in the display. The buzzer can be switched off with the **“Alarm off”** button. The **“door open”** notification appears on the display. The notification and the red light remain active until the refrigerator is closed again.

Sensor error:

If a sensor is not connected correctly or is damaged, the red LED lights up and the buzzer sounds. The buzzer can be switched off with the **„Alarm off”** button. The display shows the message **„Sensor error”**. The notification and the red light remain active until the sensor is correctly connected or replaced. The device is in emergency mode and is switched back to normal mode by acknowledgement. **Caution: Only possible when the defective sensors have been replaced or correctly connected.**

Internal temperature:

If a difference >5 degrees between the upper and lower sensor is detected or the internal temperature drops below 3 degrees, the red LED starts to light up and the buzzer sounds. The buzzer can be switched off with the **„Alarm off”** button.

The display shows the message **„Internal temperature”**. The device is in emergency mode and is switched back to normal mode by acknowledgement. **Caution: Only possible when the device is cooling in the setpoint range again.**

WARRANTY AND LIABILITY

WEPA Apothekenbedarf GmbH & Co KG (hereafter referred to as "the Manufacturer") grants the end customer (hereafter referred to as "the Customer") based a 3-year warranty from the date of delivery on the apotec® pharmaceutical refrigerator (hereafter referred to as "the Product").

The warranty covers all material and manufacturing defects. In the event of a claim made under warranty, the Manufacturer reserves the right to choose between a repair or the supply new or fully reconditioned parts. The warranty terms do not in any way affect the Customer's contractual or statutory rights.

Not covered by warranty are:

- Damage or wear to the product caused by improper or incorrect operation that is in contravention of best practice and the instructions of the Manufacturer (as laid down in this operating manual)
- Products that show signs of repairs or attempted repairs by a workshop that is not authorised by the Manufacturer to carry out such repairs
- Products that have been modified or operated with unapproved accessories
- Products whose serial number has been removed or made illegible on the type plate

To make a claim under warranty, the Customer must notify the Manufacturer prior to returning the Product by calling **+49 (0)2624 107-361** or by e-mail message to **info@wepa-apothekenbedarf.de**.

The Customer must allow the Manufacturer a period of 5 working days for the performance of a preliminary fault analysis (by telephone). Claims under this warranty shall only be considered after the Product has been returned to the Manufacturer to the following address:

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Customers based outside Germany may return the Product to their local dealer. To make a claim under warranty, please also furnish us with the original invoice bearing the date of purchase. The costs for the return of the Product shall be covered by the Manufacturer.

If a claim is made under warranty, and the inspection of the Product by the Manufacturer reveals that there is no fault, or that the damage is not covered by warranty for one of the reasons listed above, the Manufacturer shall be entitled to charge a reasonable service fee. This does not apply in cases where the Customer can prove that he could not be aware of the fact that the warranty had been voided.

The Manufacturer warranty as outlined above, including all terms and conditions, shall be transferrable to any new owner based in, provided that the original purchase invoice is submitted.

This warranty is governed by German law.

apotec® cool connect pharmaceutical refrigerator in acc. with DIN 58345

WARRANTY CARD

Warranty
 (from invoicing)
 accessories, batteries and
 consumables are precluded

Serial no.

Prod. year

Signature

Your customer no.	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	Please complete
Sender / company stamp											Date	
											Signature	

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6. Auflage: September 2021 © Copyright 2017
6th edition: September 2021 © Copyright 2017
6ème Édition : Septembre 2021 © Copyright 2017
6a edizione: Settembre 2021 © Copyright 2017
6^e editie: September 2021 © Copyright 2017
6a Edición: Septiembre 2021 © Copyright 2017