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NEDERLANDS

ESPAÑOL

apotec[®] easy pharmaceutical refrigerator in acc. with DIN 58345

USER MANUAL

apotec[®]

The professional solution for your lab.

WEPA
DIE APOTHEKENMARKE

COPYRIGHT

The copyright for this operating manual is owned by WEPA Apothekenbedarf GmbH & Co KG. This manual must not be reproduced, in any format and by any means, without the prior written permission of WEPA Apothekenbedarf GmbH & Co KG. WEPA Apothekenbedarf GmbH & Co KG shall not be liable for damage caused in connection with the use of these instructions. Errors excepted.

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TABLE OF CONTENTS

Introduction	4
Technical data	5
Safety instructions	6-11
Start-up	12-15
Troubleshooting	15-16
Warranty and liability	17
Notes	18
Warranty card	19

INTRODUCTION

The apotec® easy pharmaceutical refrigerator is used for storage of medicines requiring refrigeration and meets all the requirements of DIN 58345 such as:

- Integrated temperature recording
- Protected thermostat
- Optical and acoustic alarm
- Rechargeable battery for alarm given a power cut (12 hr. rechargeable battery)

The device meets the DIN 58345 standards only with the factory default temperature setting of 5 °C. In case of temperature adjustment (possible only by a service technician), conformity with the DIN 58345 is no longer given.

At the same time, the DIN 58345-based apotec® easy pharmaceutical refrigerator from WEPA Apothekenbedarf provides you with other benefits such as:

- 24 hours reference temperature curve at a keystroke with min./max. display
- Recirculated air cooling (see page 12)
- Voltage-free relay contact for alarm relaying
- Recirculated air cooling (see page 8)
- Automatic de-frosting and condensate evaporation (see page 11)
- Straightforward menu guide

TECHNICAL DATA

Cooling capacity	335 l
Order no.:	019861 (Door hinge on the right) 019872 (Door hinge on the left)
Power connection:	230 V / 135 W
Size:	600 x 1730 x 610 mm (W x H x D)
Weight	82 kg
Drawers:	9 drawers, max. load 15 kg per drawer (8 drawers with 2 long partitions and 6 cross-partitions, 1 drawer with 1 long partition and 4 cross-partitions)
Temperature display area:	<1°C to > 24°C Maximum and minimum temperature display
Power consumption:	0.9 kWh / 24 h

Cooling capacity	335 l – with glass door
Order no.:	019881 (Door hinge on the right) 019883 (Door hinge on the left)
Power connection:	230 V / 135 W
Size:	600 x 1730 x 610 mm (W x H x D)
Weight	89 kg
Drawers:	9 drawers, max. load 15 kg per drawer (8 drawers with 2 long partitions and 6 cross-partitions, 1 drawer with 1 long partition and 4 cross-partitions)
Temperature display area:	<1°C to > 24°C Maximum and minimum temperature display
Power consumption:	1.26 kWh / 24 h

Cooling capacity	236 l
Order no.:	019860 (Door hinge on the right) 019871 (Door hinge on the left)
Power connection:	230 V / 125 W
Size:	600 x 1323 x 610 mm (W x H x D)
Weight	65 kg
Drawers:	5 drawers, max. load 15 kg per drawer (4 drawers with 2 long partitions and 6 cross-partitions, 1 drawer with 1 long partition and 4 cross-partitions)
Temperature display area:	<1°C to > 24°C Maximum and minimum temperature display
Power consumption:	0.8 kWh / 24 h

Deliverable accessories:	Order No. 019867 Cross-partition 136 mm (sales unit: 10 pcs.) Order No. 019868 Cross-partition 206 mm (sales unit: 10 pcs.) Order No. 019869 Map cases (sales unit: 10 pcs.)
---------------------------------	--

SAFETY INSTRUCTIONS

Before using the refrigerator, read through these simple instructions !
Non-compliance might result in injury, damage to property or violation of the relevant medical safety regulations. The manufacturer shall not be liable for any damage arising from improper use of the equipment.

In general:

Check to see there is no damage to the refrigerator when unpacking it. Only operate the refrigerator as specified in the manual to ensure fail-safe operation.

Disposal pointers:

The packing is of recyclable materials.

- Corrugated cardboard/paperboard
- Mouldings of foamed polystyrene
- Polyethylene film
- Polypropylene band loops
- **The packing material is not a children's toy – risk of suffocation from the film !**
- Take the packaging to an official collection point.

Disposal

The system, except for the electronic equipment, is mainly made from steel, with some parts in aluminium. Dispose of it through recycling according to the applicable statutory waste disposal regulations. These regulations are likely to require the separate handling and disposal of

- electronic and electrical waste (printed circuit boards);
- batteries, fluorescent lamps/energy-saving bulbs, hazardous substances;
- plastics (housing);
- sheet metal, steel, copper, aluminium (separated by metal).

Before disposal, decontaminate all parts that have been in contact with process substances. Remove any hazardous materials from the the apotec® pharmaceutical refrigerator.

All oils, solutions and cleaning agents as well as contaminated cleaning tools (e.g. brushes, cloths) must be disposed of in keeping with the local requirements, in accordance with the applicable waste disposal regulations, and in compliance with the information in the material safety data sheets of the manufacturer.

Batteries:

Do not dispose of the batteries through normal household waste. You are by law obliged to ensure that batteries are processed by a specialist disposal company. To do this, bring them to the official collecting points of your municipal waste disposal service or return them to any shop selling batteries. Alternatively, you can send the batteries to us at the following address:

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We will dispose of them properly for you without any charge. Batteries containing toxic substances are marked with a sign consisting of a crossed out dustbin and the chemical symbol (Cd, Hg or Pb) indicating the hazardous substance (heavy metal). The crossed-out waste bin symbol indicates that rechargeable and non-rechargeable batteries must not be emptied into the normal household waste. The chemical symbols indicate the following: Pb: battery contains lead; Cd: battery contains cadmium; Hg: battery contains mercury.

Place of installation / Power connection:

- Protect the refrigerator from direct sunlight!
- The floor must be level and horizontal at the location. Offset any unevenness with the adjusting feet.
- Ensure there is adequate ventilation and venting!
- Do not use an extension lead to connect the refrigerator with other appliances – risk of over-heating!
- The place of installation must have a temperature range from 10 °C to 35 °C. The relative humidity may not exceed 75%.

Power connection:

Type of current (AC) and voltage at the point of installation must be the same as the particulars on the type plate. The type plate is inside the refrigerator on the left-hand side. Fuse the socket at 10 amperes or higher. It must be away from the appliance back panel and be easy to access. Only connect the refrigerator by way of a properly earthed earthing-contact socket.


Recirculated air cooling:

Recirculated air cooling involves the swirling of cooled air with a fan. This process enables the cold air to be rapidly and evenly spread inside. Hardly any temperature stratification occurs; the temperature is practically the same throughout the entire interior.

Safety and warning instructions:

- To avoid any injuries or material damage, two people should be entrusted with unpacking and setting up the refrigerator.
- If any damage is noticed on the refrigerator, immediately contact the supplier – before connecting the device.
- To ensure fail-safe operations, only install and connect the refrigerator in accordance with the instructions in the manual.
- If a fault occurs, disconnect the refrigerator from the mains, pull out the mains plug or trip/remove the fuse. **To disconnect the refrigerator from the mains, pull the mains plug itself and not the connecting cable.**
- **Interventions and repairs to the refrigerator must be carried out** by the competent after-sales service. Failure to do so creates considerable risks for the user. The same applies to the replacement of the power connection line.
- **Inside the refrigerator**, do not undertake work involving any naked flame or sources of ignition.
- Ensure that the refrigeration circuit is not damaged when handling or cleaning the refrigerator. Keep sources of ignition away from any damaged points and ventilate the area well.
- **Do not use base, drawers, doors etc.** for stepping on or for propping up.
- This refrigerator is not intended for use by people (including children) with physical, sensory or mental impairments or for those who are inexperienced in its use or who lack the necessary knowledge – unless they have been instructed in using the refrigerator by someone who is responsible for their safety or who supervises them in the initial stages. Children should be supervised at all times to ensure they do not play with the refrigerator.
- Avoid constant skin contact with cold surfaces or cooled/frozen products. This can otherwise lead to pain, numbness and frost-bite. Provide for protection if work involves lengthy periods of skin contact e.g. by wearing gloves.

- Do not store in the refrigerator any explosive materials or spray cans with combustible propellants such as propane, butane and pentane. Any escaping gases could be ignited by the electrical components. These kinds of spray cans can be recognized as such by the printed details on content or a flame symbol
- Do not use any electrical equipment inside the refrigerator.
- Keep keys of lockable equipment well out of the reach of children.
- Only load the refrigerator with 2 - 8 °C refrigerated products (intrinsic temperature). Proper cooling cannot be ensured if this is disregarded!
- Do not put any foodstuffs/drinks into the refrigerator!
- Visually check information labels, type plate and magnetic contact at least twice a year as to their legibility/intactness. If any damage is detected, please immediately contact the after-sales service.
- Regularly check (including on weekends) the refrigerators as to their working order. Regularly means every 24 hours. (Recommendation GSM module; errors are reported immediately to one or more mobile phones in case of failure – order No. 019877).
- Do not keep any "frozen products" (cold packs/pads, ice etc.) in the refrigerator. Proper cooling cannot be ensured if this is disregarded!
- To ensure a proper data record, date and time must always be up to date! Please note summer and winter time.
- The internal memory records the operating data for 10 months and is then overwritten. Please back up the data on a regular basis.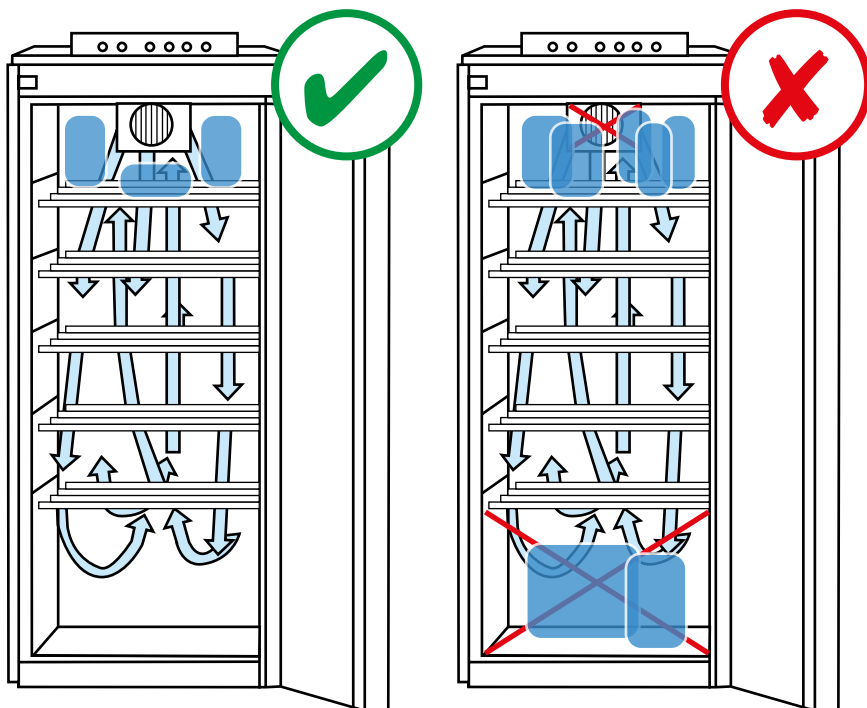

The lower/upper section of the refrigerator (see drawing) must not be blocked. The recirculation function will be disturbed and ventilation cannot be guaranteed!

- The mains plug must always be freely accessible!
- The device is not suitable for installation, e.g. in a cabinet.
- The electronic unit with display must be easily accessible, to be able to operate the rear panel if necessary.

WARNING:

Risk of injury due to electric shock! The device may only be opened by the customer service or specially trained personnel.

- According to the EN 378 standard, the room where your device is installed must have a volume of 1 m³ per 8 g of refrigerant R 600a, so that no ignitable gas-air mixture can form in the installation room in the event of a leak in the refrigerant circuit.
The refrigerant quantity is indicated on the rating plate inside the device.

CAUTION:

The device must be aligned horizontally and vertically. If the device is tilted, the body can become deformed and the door will not close properly. Level out uneven floors with the adjustable feet.

- Do not connect the device via extension cables or distribution sockets.

Fault:

You can correct the following faults yourself by checking the possible causes:

- **The appliance does not work:** Please check
 - if the refrigerator is switched on,
 - if the mains plug fits correctly in the socket
 - if the socket fuse is OK
- **The noise emission is too loud:** Please, check if
 - the refrigerator rests firmly on the floor,
 - the operating refrigerating unit is the reason for vibrations in furniture or objects next to it.Note that flow noises in the refrigeration circuit cannot be prevented.

If none of the above causes apply and you cannot correct the fault yourself, then please contact the after-sales service.

De-frosting:

The cooling section defrosts automatically.

If the rear wall still freezes over, the unit must be defrosted manually.

- To defrost switch off the refrigerator. Disconnect the mains plug
- Remove the cooling products from the refrigerator and store where cool.
- Leave the refrigerator door open when defrosting. Take up any condensed water with a cloth and clean the refrigerator.

When defrosting do not use any mechanical implements or other artificial aids save for those recommended by the manufacturer.

Cleaning:

- Place the refrigerator out of service before cleaning (pull out the mains plug).
- Clean the interior, fittings (drawers) and external panels with tepid water and a little rinsing agent. Do not use any cleaning or chemical solvents containing sand or acid.
- Do not work with steam cleaning devices Risk of damage/injury
- Make sure that no cleaning water gets into the electrical parts or into the ventilation grate.
- The chiller with the heat exchanger (metal grate at the back of the refrigerator) should be cleaned or dusted once a year.
- Clean the drain in the refrigerator with something thin (Q-tips and the like)

OUTER CONTROL PANEL BUTTONS – EXPLANATION OF TERMS, COMMISSIONING AND FUNCTIONS

Menu button:

call up the menu or if you are in the menu, go one step back in the menu.

OK button:

Button for acknowledgement in the menu.


Arrow Up button:


navigate upward (scroll) in the menu.

Arrow downbutton: scroll down in the menu.

SD card slot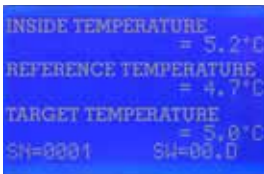

The refrigerator provides an easy way to display the current temperature at a glance. To do so, press "OK" while the status screen is active.

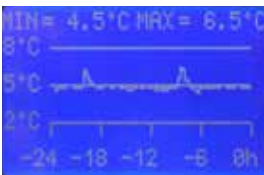In addition, you can view reference temperature gradient and the min./max. reference temperatures of the last 24 hours. While the status screen is active, press arrow "up" or "down".

Note:

You will notice a short-term temperature increase every day, which is due to the automatic defrosting process. Here the 8 °C reference temperature is usually not exceeded in the normal case.

Internal temperature:

Actual measured air temperature in the refrigerator.

Reference temperature:

"Simulates" the temperature of the refrigerated products. In accordance with DIN 58345, this is recorded in a built-in reference body and appropriately adjusted. This temperature is also decisive for the alarm.

Target temperature:

The target temperature is preset in accordance with DIN 58345 at 5 °C. A deviation of ± 3 °C is permitted. The deviation refers to the reference temperature. If the temperature range is exceeded or undershot, the refrigerator emits an optical and acoustic alarm.

LEDs:

In addition to the indicated errors or alerts in the display, the LED lamps show an overview of the operating status:

Green:	Proper refrigerator operation
Yellow:	Battery weak, data backup necessary, menu mode, short power failure (under 5 seconds)
Red:	Power failure, temperature error (reference temperature), sensor defective

**Start-up:**

The refrigerator is in the non-operating mode at the time of delivery: The display is non-active. As soon as you connect the mains plug, you turn off the sleep mode. A rechargeable battery integrated in the refrigerator needs to be charged for 36 hours on first commissioning.

Menu:

You can make the following settings in the menu.

- Read SD card
- Set time
- Set date
- Switch off
- Set language

In addition, there is a service menu that is secured by a PIN and is only accessible by the customer service!

Setting the time:

To set the time, start by pressing the "Menu" button. Now go to the "Time" menu item "Set Time". Push the button "OK". To set the hour, scroll with the arrow keys to the desired time (hour). Now repeat the process to set the minutes. When you have set the minutes, please confirm with "OK". The menu will be automatically exited.

Setting the date:


Start by pressing "Menu" to access the menu. Now use the arrow keys to change to the menu item "Set Date" Press "OK" to continue. Now proceed as with setting the time. First the year, then the month and finally the day. Confirm the day by pressing "OK". The menu will be automatically exited.

Data storage on SD card:

The internal temperature and reference temperature are stored on the internal memory of the device every 60 seconds. In addition, special events are stored, such as fluctuations in temperature, reference temperatures, acknowledgements, etc. as well as the opening and closing of the door. The apotec® pharmaceutical refrigerator provides a function to transfer this data from the memory to an SD card and read it on a PC with the supplied apotec® refrigerator documentation. For this, proceed as follows:

Insert your SD card into the envisaged slot. Now activate the "Menu" and scroll to the point "Data backup". Press "OK" to transfer the information. The percentage progress will be displayed.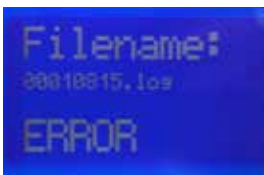

If now no SD card is in the slot, "Error" appears on the display and the menu will be automatically exited. Depending on data volume, the operation takes up to 10 minutes. After the transfer, the SD card can be removed and data can be read on the PC.

Setting the language:

To set the language, select the menu item "Language" – press "OK". Now please Select the desired language and confirm with "OK".

Switching off:

You can turn the unit off manually. To do this, select the menu item "Off" and confirm with "OK".

TROUBLESHOOTING

Power failure:

If the refrigerator is disconnected from the mains cable or no power gets to it (due to a power cut), then this fault message "Power failure" and the **backlight is dimmed**. (Note: The display remains in this state for min. 12 hours until power supply is available again). The green lamp signalling smooth refrigerator function goes out. Subsequently, the red LED lights up and the built-in buzzer sounds. In this case, the buzzer can be disabled with the button "OK". Once the power supply or line voltage have been re-established, the error message disappears after acknowledgement with "OK" and normal operation is continued.

Reference temperature limit values:

If the factory set temperature of 5 °C is exceeded or undershot by 3 °C, the red LED lights up and the built-in buzzer sounds. The buzzer can be switched off by using the "OK" button. The message "Reference temperature" appears on the display. The message and the red lamp remain active until the refrigerator is again within the target range.

Door open:

If the door is open longer than 120 seconds, an optical and acoustic alarm will be triggered. You can disable the buzzer by pressing the "OK" button. The message "Door" will appear on the display. The message and the red lamp remain active until the refrigerator is closed again.

BATTERY:

If the battery does not have sufficient capacity, the error message "AKKU" appears.

Note: Since the rechargeable battery is a wearing part, it is not covered by the warranty/guarantee.

The battery should then be replaced by an authorised service centre.

Otherwise the alarm function and data recording for at least 12 hours cannot be guaranteed in case of power failure.

Sensor error:

If a sensor is not connected correctly or is damaged, the red LED lights up and the buzzer sounds. The display shows the message "Sensor error".

The notification and the red light remain active until the sensor is correctly connected or replaced. The audible alarm can be suppressed with the "OK" key.

WARRANTY AND LIABILITY

WEPA Apothekenbedarf GmbH & Co KG (hereafter referred to as "the Manufacturer") grants the end customer (hereafter referred to as "the Customer") based a 3-year warranty from the date of delivery on the apotec® pharmaceutical refrigerator (hereafter referred to as "the Product").

The warranty covers all material and manufacturing defects. In the event of a claim made under warranty, the Manufacturer reserves the right to choose between a repair or the supply new or fully reconditioned parts. The warranty terms do not in any way affect the Customer's contractual or statutory rights.

Not covered by warranty are:

- Damage or wear to the product caused by improper or incorrect operation that is in contravention of best practice and the instructions of the Manufacturer (as laid down in this operating manual)
- Products that show signs of repairs or attempted repairs by a workshop that is not authorised by the Manufacturer to carry out such repairs
- Products that have been modified or operated with unapproved accessories
- Products whose serial number has been removed or made illegible on the type plate

To make a claim under warranty, the Customer must notify the Manufacturer prior to returning the Product by calling **+49 (0)2624 107-361** or by e-mail message to **info@wepa-apothekenbedarf.de**.

The Customer must allow the Manufacturer a period of 5 working days for the performance of a preliminary fault analysis (by telephone). Claims under this warranty shall only be considered after the Product has been returned to the Manufacturer to the following address:

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Customers based outside Germany may return the Product to their local dealer. To make a claim under warranty, please also furnish us with the original invoice bearing the date of purchase. The costs for the return of the Product shall be covered by the Manufacturer.

If a claim is made under warranty, and the inspection of the Product by the Manufacturer reveals that there is no fault, or that the damage is not covered by warranty for one of the reasons listed above, the Manufacturer shall be entitled to charge a reasonable service fee. This does not apply in cases where the Customer can prove that he could not be aware of the fact that the warranty had been voided.

The Manufacturer warranty as outlined above, including all terms and conditions, shall be transferrable to any new owner based in, provided that the original purchase invoice is submitted.

This warranty is governed by German law.

apotec® easy pharmaceutical refrigerator in acc. with DIN 58345
WARRANTY CARD

Warranty
 (from issuing the invoice)
 Accessories and consumables excluded!

Serial No.

Prod. Year

Signature

Your customer no.	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	Please complete
Sender / company stamp											Date	
											Signature	

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5. Auflage: September 2021 © Copyright 2015
5th edition: September 2021 © Copyright 2015
5e Édition : Septembre 2021 © Copyright 2015
5a edizione: Settembre 2021 © Copyright 2015
5^e editie: september 2021 © Copyright 2015
5ta Edición: Septiembre 2021 © Copyright 2015